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21st century is the era of biotechnology. Biotechnology has been becoming a vital approach to solve the emerging problems of human subsistence. Biotechnology now is playing a more and more important role in varied fields, such as bioenergy, biomaterial, environment protection, biopharmaceutical, modern agriculture, and modern food, etc.

High-efficient bioreactors (fermenters) are essential for the industrialization of biotechnology. Eastbio is a professional company devoted to the industrialization of biotechnology.

Eastbio focus itself on the whole production process of bio-products, from raw material treatment, core bioreactor, down process and bioprocess automatization, to wastewater treatment of the fermentation plant. Our belief is that the equipment should satisfy the requirements of process and product.

Our vision is "Be efficient and honest, be science and technology-oriented". Our base is "Be specialized and focused, provide technology and service".

In 1996, Eastbio was established by the professional researchers from Jiangsu University. And today Eastbio has become the most famous provider of professional equipment for biotechnology industrialization in China.

We hope to become the friend of our customers and grow with you as well as create value for you.

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Ninth CPPCC National Committee of Jiangsu Province

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Ninth CPPCC National Committee of Zhenjiang City

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Visiting Professor of Tianjin University Science and Technology

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Enzyme Engineering Professional Committee of China

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Editorial Board Member of Journal of China Pharmaceutical University

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Zhenjiang East Biotech Equipment and Technology Co., Ltd. which was established in 1996, adopted "highly efficient, integrity & technology-based" as business philosophy, persists in servicing the field of biological engineering as to create extra value for customers.

Eastbio has always focused in bioengineering area and commits itself to industrialize biotechnology, as well as providing total bioengineering solution for customers. Eastbio has grown and becoming the most famous manufacturer of biotechnological equipments in China until today. The trademark of "Eastbio" has also become a byword for quality and excellent in terms of product and service.

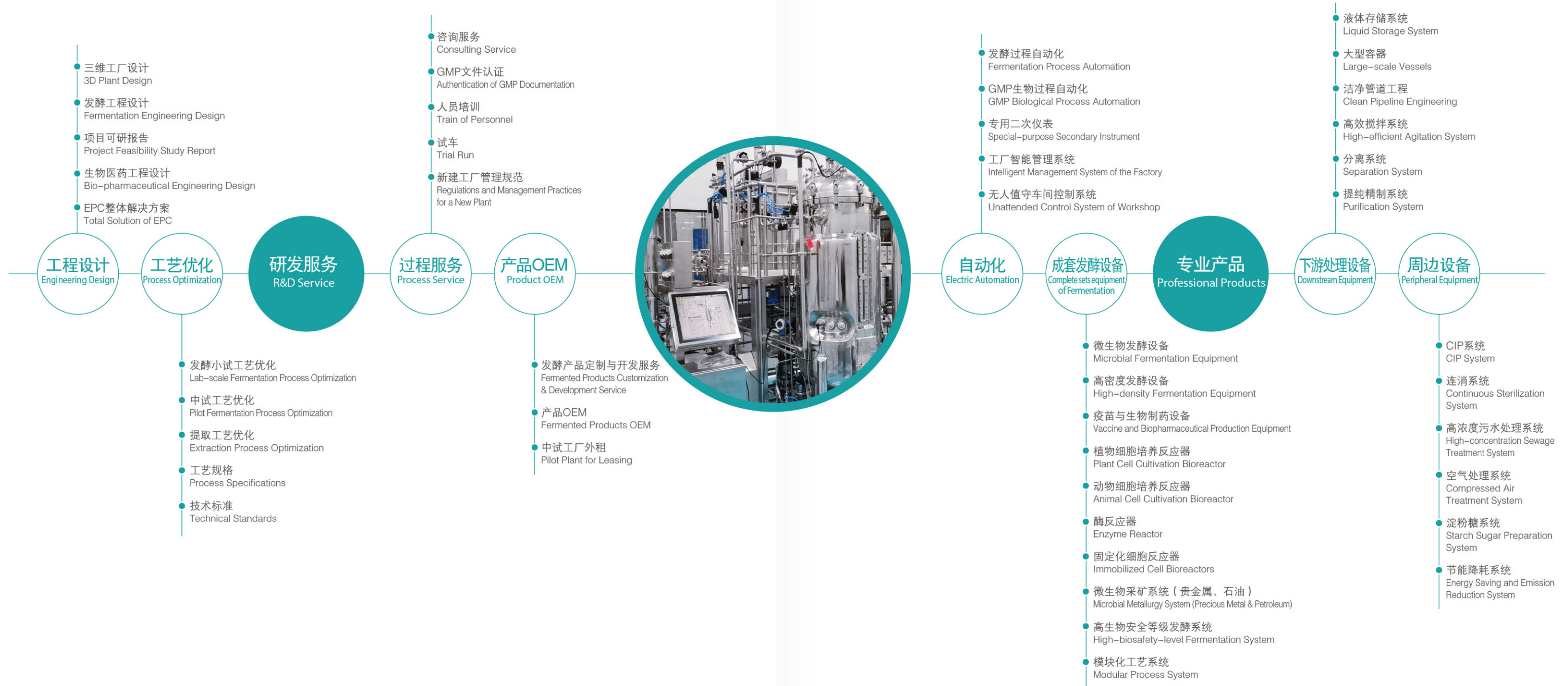
d) Besides, the company is currently cover a total area of 153,000 square meters, including the plant area of 100,000 square meters, with a total registered capital of RMB 60 million Yuan, has capabilities in R & D, design, manufacturing and integrated services.

Jiangsu Multifunctional Fermentation Equipment Engineering Research Center (JMFERC) and Zhenjiang Bioengineering Research Center (ZBRC) are affiliates of the company.

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Fermentation Project Turn-key Contractor

型号(Model)	GBJSXN/GBCSXN / 3 / 5 / 8
公称体积(Nominal Volume)	3L / 5L / 8L

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#### Application Scope:

It is an ideal tool for precise fermentation experiment in microbiological laboratory for both research centers, institutes as well as private enterprises. It is normally applied for culture medium selection, fermentation process optimization and verification of strains as well as its production process.

#### Technical Specification:

Loading Quotient	65-80%
Material	Special autoclavable glass
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Magnetic stirred from bottom / Mechanical stirred from top end
Sealing Form	iii.Magnetic coupling without mechanical seal / Mechanical seal
Structure	Cabinet type / Benchtop
Sterilization	Off-situ sterilized in an autoclave
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. It applies magnetic coupling to eliminate mechanical seal leakage risk. The fermenter can be observed clearly, dismantled and mounted easily, and operated conveniently.

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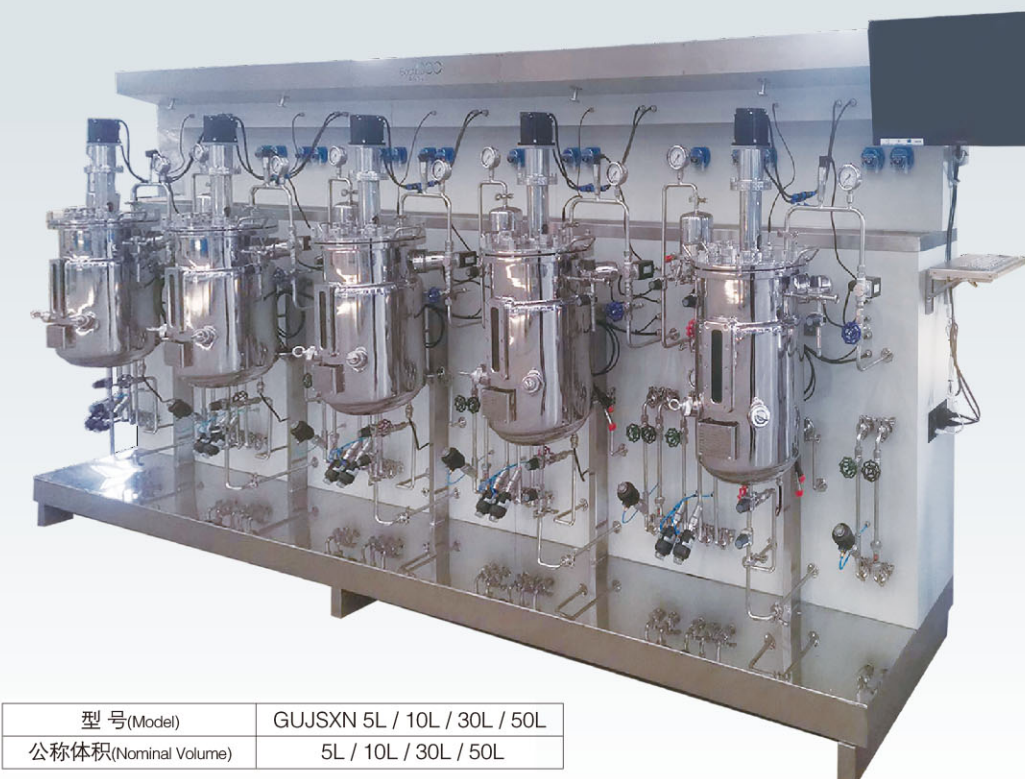
Technical Specification:

Loading Quotient	65-80%
Material	Special autoclavable glass
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Mechanical stirred
Sealing Form	Special mechanical seal with reliable performance
Structure	Benchtop
Sterilization	Off-situ sterilized in an autoclave
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	It is more suitable for sifting through culture medium, optimizing process and comparing properties of strain as multiple vessels run simultaneously to perform parallel test. The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. High quality special mechanical seal is applied to assure good gas tightness. The inside situations can be observed clearly, and the assembled cabinet are able to put in use conveniently.

型号(Model)	GUJSXN 5L / 10L / 30L / 50L
公称体积(Nominal Volume)	5L / 10L / 30L / 50L

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Technical Specification:

Loading Quotient	65-75%
Material	SUS316L/SUS304 (glass shell can be applied to fermenters with volume below 10 liters)
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Magnetic stirred from bottom
Sealing Form	Special mechanical seal with reliable performance
Structure	Floor type with European style frame
Sterilization	In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The in-situ sterilization process is similar to the process of scale-up experiments. The operation method is simple and the sterilization is safe and reliable. Special mechanical seal is applied to assure good gas tightness. The inside situations can be observed clearly via wide viewing glass. The equipment can be dismounted and mounted easily. Frameset construction is convenient for operation.

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
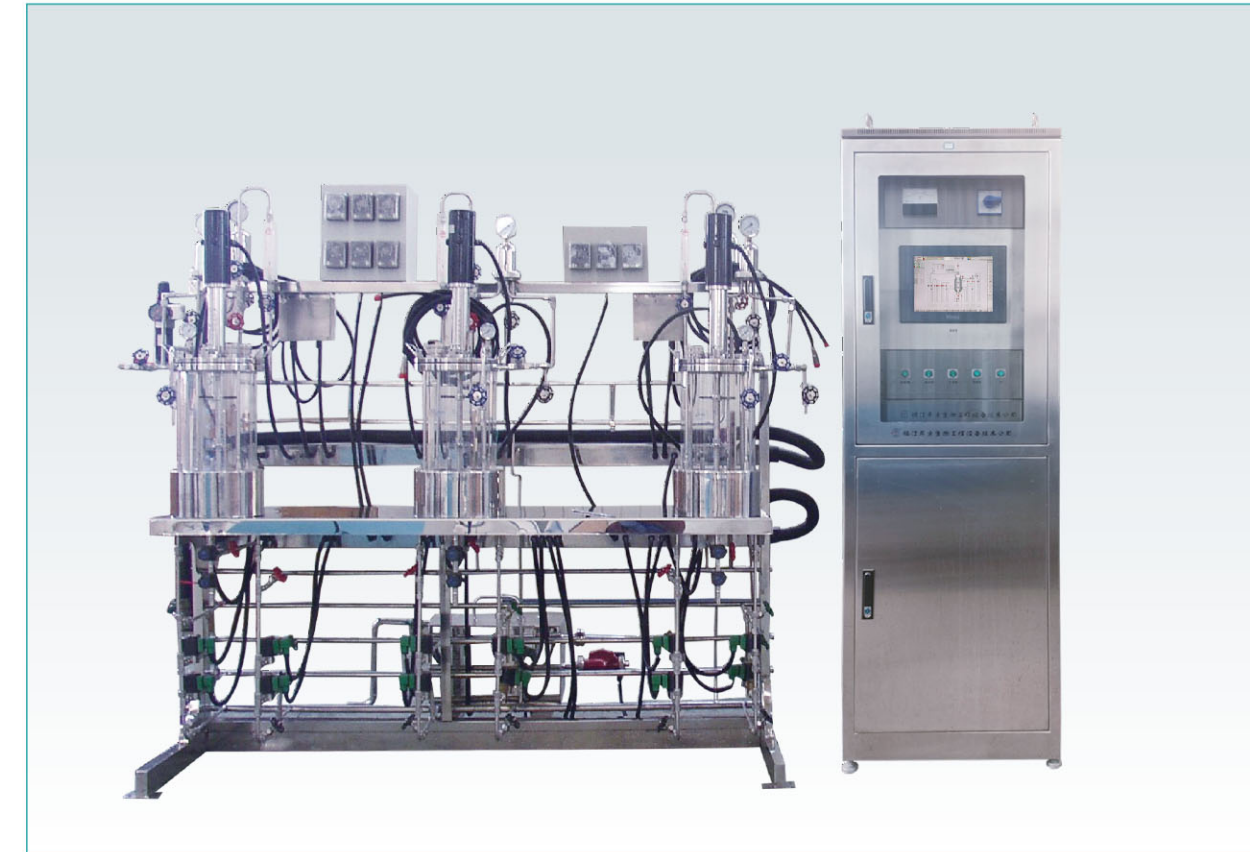
#### Application Scope:

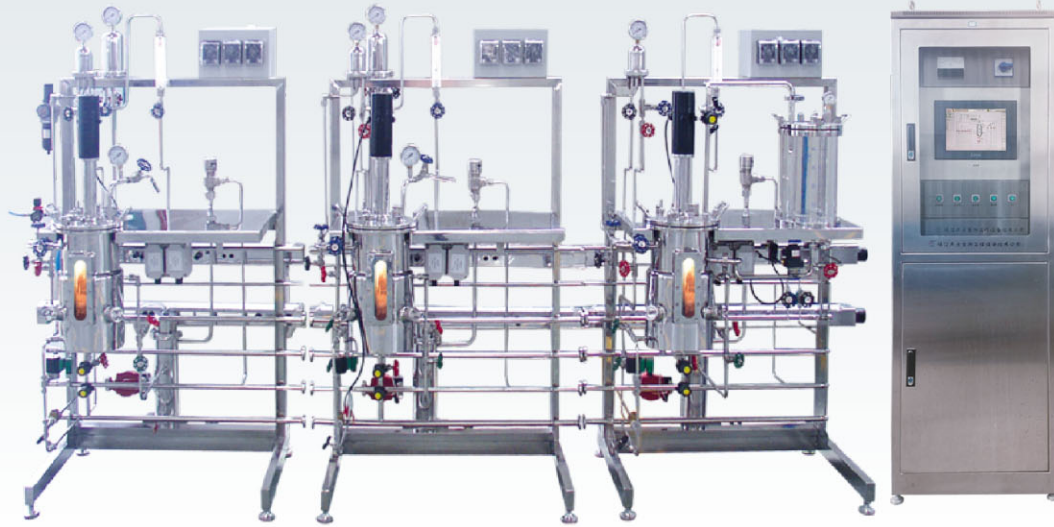It is an ideal tool for precise fermentation experiment in microbiological laboratory for both research centers, institutes as well as private enterprises. It is normally applied for culture medium selection, fermentation process optimization and verification of strains as well as its production process. The incorporated in-situ sterilization technology has also ease the up scaling due to comparable fermentation result.

#### Technical Specification:

Loading Quotiety	65-75%
Material	Special autoclavable glass
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	ii.Magnetic stirred from bottom / Mechanical stirred from top end
Sealing Form	Magnetic coupling without mechanical seal / Mechanical seal
Structure	Floor type
Sterilization	v.In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. It applies magnetic coupling to eliminate mechanical seal leakage risk. The fermenter can be observed clearly, dismantled and mounted easily, and operated conveniently.

型号(Model)	GUJS 5L / 50L / 100L
公称体积(Nominal Volume)	5L / 50L / 100L

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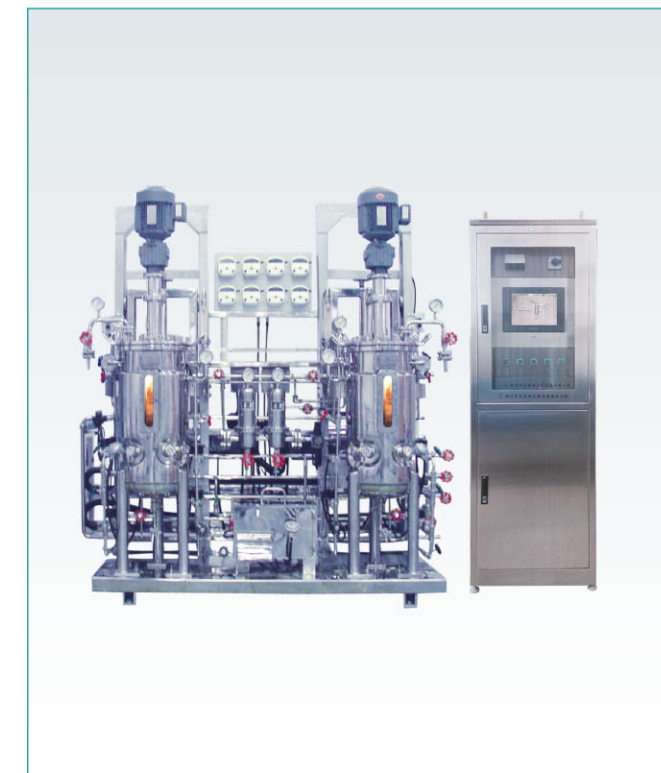
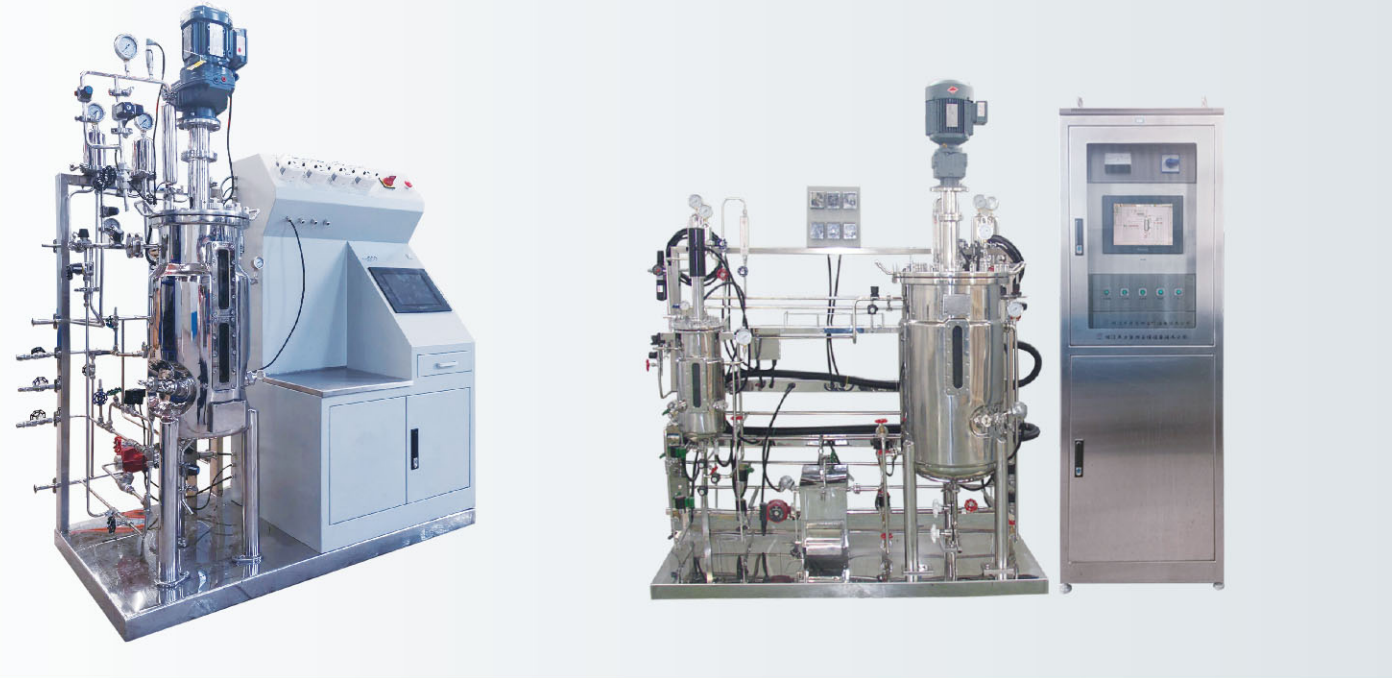
#### Application Scope:

It is an ideal tool for precise fermentation experiment in microbiological laboratory for both research centers, institutes as well as private enterprises. It is normally applied for culture medium selection, fermentation process optimization and verification of strains as well as its production process. The incorporated in-situ sterilization technology has also ease the up scaling due to comparable fermentation result.

#### Technical Specification:

Loading Quotient	65-75%
Material	SUS316L/SUS304
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Mechanical stirred from top end
Sealing Form	Mechanical seal
Structure	Floor type
Sterilization	In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. It applies magnetic coupling to eliminate mechanical seal leakage risk. The fermenter can be observed clearly, dismantled and mounted easily, and operated conveniently.

型号(Model)	GUJT / 100 / 300 / 500
公称体积(Nominal Volume)	100L / 300L / 500L

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#### Application Scope:

It is an ideal tool for precise fermentation experiment in microbiological laboratory for both research centers, institutes as well as private enterprises. It is normally applied for culture medium selection, fermentation process optimization and verification of strains as well as its production process.

#### Technical Specification:

Loading Quotient	65-80%
Material	SUS316L/SUS304
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Mechanical stirred from top end
Sealing Form	Mechanical seal
Structure	Floor type
Sterilization	In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. It applies magnetic coupling to eliminate mechanical seal leakage risk. The fermenter can be observed clearly, dismantled and mounted easily, and operated conveniently.

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**Application Scope:**

It is an ideal tool for precise fermentation experiment in microbiological laboratory for both research centers, institutes as well as private enterprises. It is normally applied for culture medium selection, fermentation process optimization and verification of strains as well as its production process.

**Technical Specification:**

Loading Quotient	65-80%
Material	SUS316L/SUS304
Agitation	Flat-blade/oblique blade/curved blade/axial flow impeller for choice
Type of Drive	Mechanical stirred from top end / Magnetic stirred from bottom
Sealing Form	Mechanical seal / Magnetic coupling without mechanical seal
Structure	Floor type
Sterilization	In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. It applies magnetic coupling to eliminate mechanical seal leakage risk. The fermenter can be observed clearly, dismantled and mounted easily, and operated conveniently.

型号(Model)	Pr/GUE / 10L / 1000L / 10000L
公称体积(Nominal Volume)	10L / 1000L / 10000L

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#### Technical Specification:

Loading Quotient	65-80%
Material	Borosilicate glass / SUS316L/SUS304
Type of Drive	Airlift type
Structure	Floor type
Sterilization	In-situ sterilization
Control method	DF type control system based Rigour™ integrated controller, displayed and operated on touch screen.
Characteristics	It is more suitable for sifting through culture medium, optimizing process and comparing properties of strain as multiple vessels run simultaneously to perform parallel test. The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. High quality special mechanical seal is applied to assure good gas tightness. The inside situations can be observed clearly, and the assembled cabinet are able to put in use conveniently.

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Application Scope:

It is widely used in microbiological laboratory of scientific research institutes and enterprises for solid state fermentation experiments. And it is best choice for solid state fermentation experiments and small scale production, especially suit solid state fermentation of raw materials of medicines and health products.

型号(Model)	SF / 50 / 100
公称体积(Nominal Volume)	50L / 100L

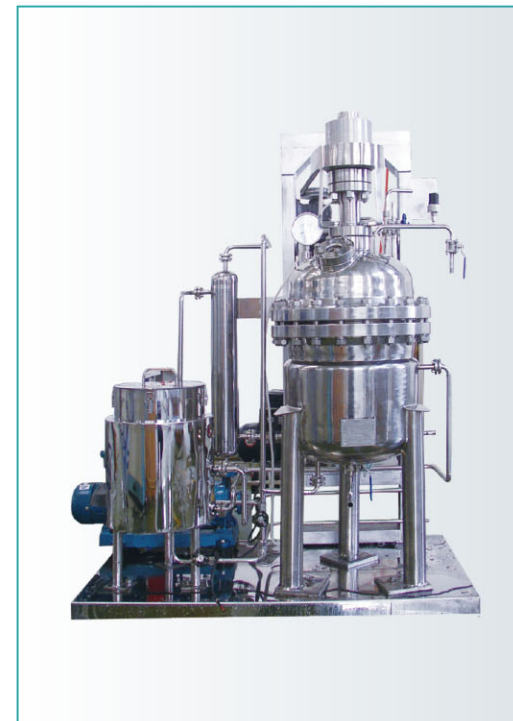
型号(Model)	GUM
公称体积(Nominal Volume)	7L / 10L / 50L / 500L.....50KL

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Technical Specification:

Loading Quotient	20-55%
Material	Stainless steel
Agitation	rotary-shaft stirrer
Type of Drive	Mechanical stirred from side part
Sealing Form	Special mechanical seal with reliable performance
Sterilization	In-situ sterilization
Control method	Instrument/computer control. The system can automatic control and real-time display parameters as temperature, humidity, pressure, flow rate, etc., and can record, analyze and print the data too.

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Application Scope:

It is ideal tool for scientific research institutes or enterprises in mineral study and production. It can be applied to rare metal, precious metal or nuclear reactive metal enrichment purpose as well as handling of environment by metal contamination. The size ranges from laboratory to production scale.

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Technical Specification:

Loading Quotient	65-70%
Material	SUS316L
Agitation	Special impellers or combination impellers
Type of Drive	Mechanical stirred from top. Self-balancing device is provided to the large reactor.
Sealing Form	Special mechanical seal with reliable performance
Sterilization	In-situ sterilization
Control method	Ebio™ control system based SIEMENS PLC, displayed and operated on touch screen.
Characteristics	The equipments match the process perfectly through special design according to certain process. The sterilizing operation is safe and reliable, and the substrate concentration is accurate after sterilization. The inside situations can be observed clearly via wide viewing glass. The equipment can be dismantled and mounted easily. Special mechanical seal is applied to assure good gas tightness. The agitation system is powerful for adequately mixing.

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Application Scope:

It is ideal tool for plant cell cultivation and widely used in microbiological laboratory for scientific research institutes and enterprises. It can be applied to plant cell cultivation and metabolite production for yew cell, vinca cell, tobacco cell and ginseng cell. It ranges from laboratory to production scale.

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Technical Specification:

Loading Quotient	65-80%
Material	SUS316L (glass shell can be applied to 5L fermenter)
Agitation	p-mix type special impellers
Type of Drive	Mechanical stirred from top with rapid dismounting stirrer
Sealing Form	Special mechanical seal with reliable performance
Structure	Frame style
Sterilization	In-situ sterilization (5L can be autoclaved)
Control method	Ebio™ control system based SIEMENS PLC, displayed and operated on touch screen.
Characteristics	The sterilizing operation is safe and reliable, and the culture medium concentration is accurate after sterilization. Multiple nutrient solutions supply. Special agitation system is applied to satisfy biological need of plant cell. The inside situations can be observed clearly via wide viewing glass. It can be dismounted and mounted easily and operated conveniently.

型号(Model)	CAJS / 5 / 10 / 50 / 100
公称体积(Nominal Volume)	5L / 10L / 5L / 100L

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Application Scope:

It is generally used in scientific research institutes and enterprises for cell suspension culture as well as microcarrier adherent culture. It facilitate scale-up of the cultivation technology, and it can provide optimal conditions to the growth and proliferation of animal cell. It ranges from laboratory to production scale.

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Technical Specification:

Loading Quotient	70-80%
Material	SS316L (glass vessel can be applied to 5L & 10L bioreactor)
Agitation	Imported special bottom magnetic agitation system, particular impellers for mammalian cell cultivation
Type of Drive	Four-line gases mixing & permeating system
Control method	V2011 control system with Siemens PLC and HMI (Touch screen monitor)

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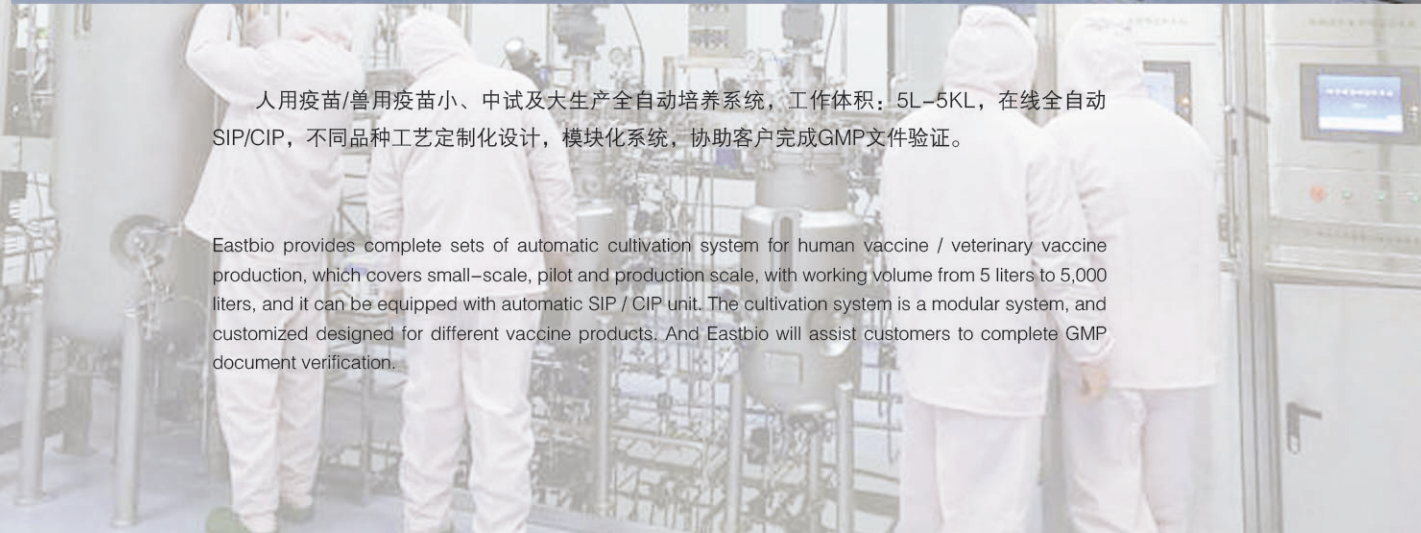
Application Scope:

It is widely used for pharmaceutical industry in scientific research and industry production. Eastbio is capable to be contractor on turnkey basis including design, fabrication, installation, commissioning, training and technological system. In addition, Eastbio could assist customer in carrying out various kinds of GMP certification based on customer requirement.

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Eastbio provides complete sets of automatic cultivation system for human vaccine / veterinary vaccine production, which covers small-scale, pilot and production scale, with working volume from 5 liters to 5,000 liters, and it can be equipped with automatic SIP / CIP unit. The cultivation system is a modular system, and customized designed for different vaccine products. And Eastbio will assist customers to complete GMP document verification.

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Application Scope:

It's the overall solution for vaccine preparation system. Eastbio can provide engineering general contracting including process design, fermentation module, buffer / medium preparation module, purification module, harvest module, CIP module and other auxiliary process equipment, and assist users in various GMP qualification demonstrations.

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Application Scope:

It is widely used in the production of probiotics such as Bifidobacterium / Enterococcus faecalis / Lactobacillus. Eastbio can provide engineering general contracting including design, manufacturing, installation, commissioning, personnel training, process schedule, etc., and can assist users to conduct various GMP qualification demonstration in combination with target products.

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Application Scope:

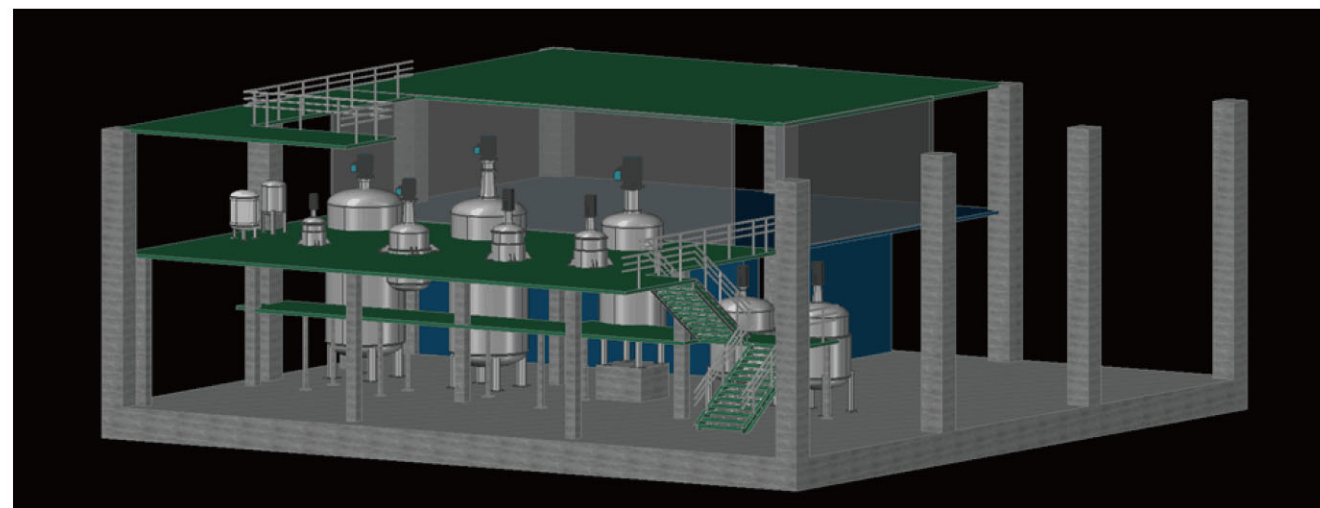
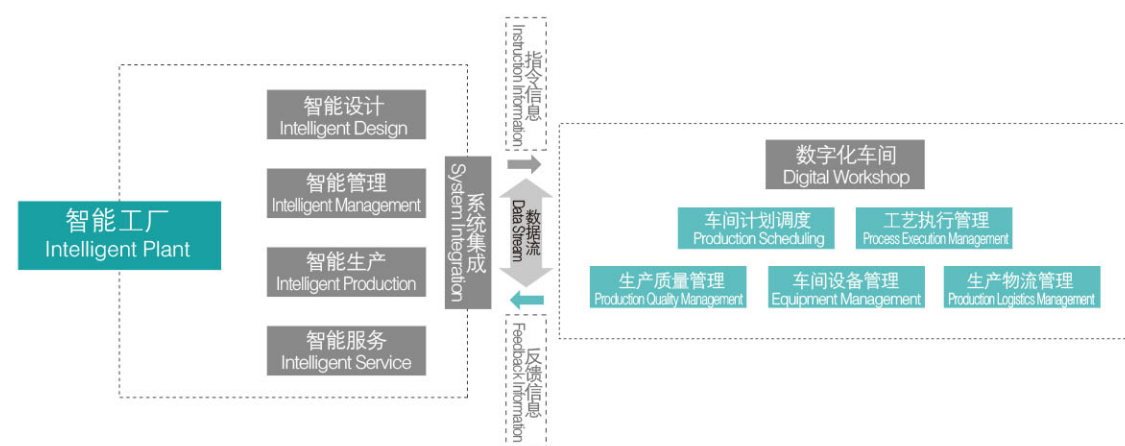
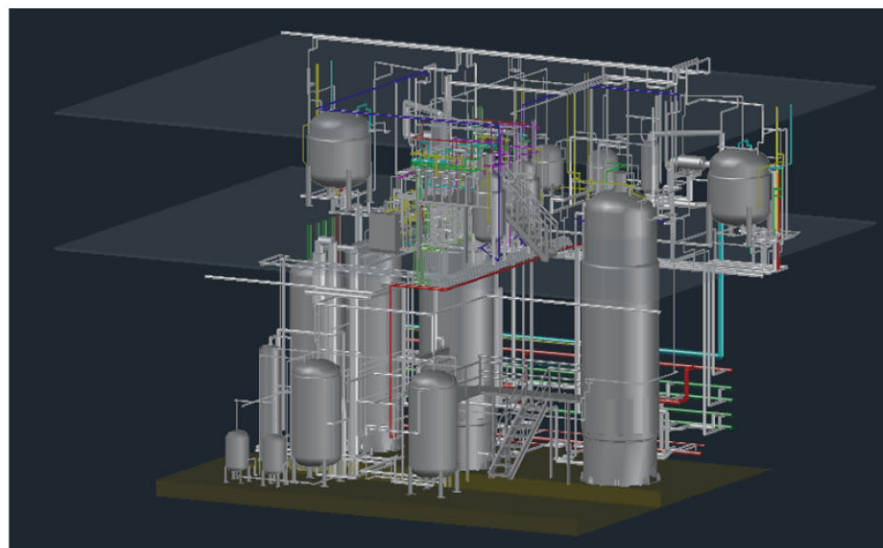
It is suitable for automatic intelligent factory and GMP biopharmaceutical project.

The bottom control layer (fermentation system, material process, etc.) is the field unit module components, which adopts PLC / DCS automatic control management unit.

The middle level is the factory equipment intelligent management software (MES), data link up and down transmission equipment operation automatic control software, equipment operation management software, equipment automatic maintenance management software, etc.

The upper plan layer is the factory data management system, which adopts ERP system to detect and feedback the production, supply and sales cost management of the whole plant, including financial management, production management, sales management, comprehensive management, operations management, etc.

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Eastbio systematically design the complete fermentation plant with the international advanced intelligent 3D plant design system, which covers:

- Intelligent PID
- Equipment module
- Steel structure
- Three dimensional pipeline
- Utilities
- Ventilation pipes HVAC
- Cable bridge
- Electric automation
- Continuity checking
- Collision checking
- 3D model roaming checking……

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Zhenjiang East Biotech Equipment and Technology Co., Ltd. has well qualified manpower and comprehensive capability to undertake fermentation engineering turnkey project. Eastbio has vast experience in microbial fermentation technology research as well as scaling-up, technological advanced design capacity, advanced machine manufacturing ability, integrated project organization and execution, commissioning and personnel training capability. Eastbio has gained a good reputation & recognition in bio-processing market by executing several fermentation engineering turnkey projects in Asia successfully.

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Eastbio adopts the world-wide common used DCS system to develop the automatic control system of biological fermentation process and biological fermentation plant, which can analyze and control up to 13 parameters.

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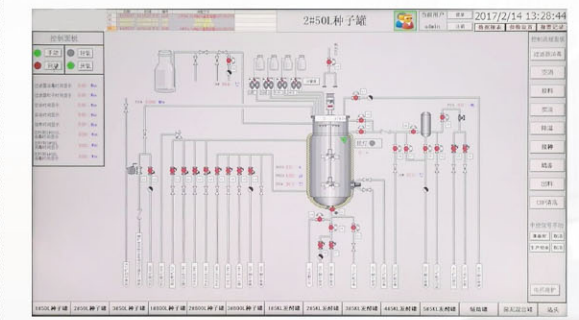
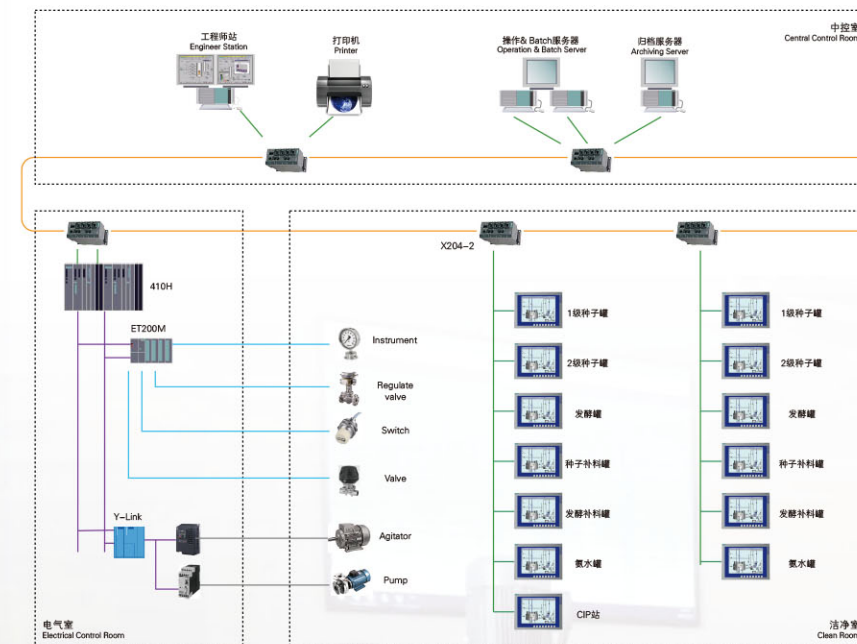
EBIO™ industrial biochemical process automatic control system is a new type of control system that base on industrial scale PLC which was developed by our in-house and external experts for a duration of 3 years. As it incorporates industrial features, it can control each parameter in the fermentation process that meeting physiological need for the best cell growth. It controls up to 12 fermenters at a time with single controller.

The system can perform set-point control, sequence control and program control. And it can also perform various kinds of comprehensive data analysis, auto-reports generation and results analysis. The interface is visually, and the touch screen can display process flow chart, working conditions for each parameter. The system can realize automatic Sterilization-in-Place (SIP) too.

There are 13 parameters for customers to choose.

The system can be extended for the entire plant automatic control including media preparation, fermentation, extraction, purification and warehouse management.


The parameters that can be measured and controlled include temperature, agitation speed, foam, pH, DO, pressure, air flow rate and multiple routes feedings, and we can also test 12 optional parameters such as oxidation-reduction potential, glucose, lactic acid and special amino acid according to the requirements of customers. Customers can select the optional parameters based on their respective experimental requirement.

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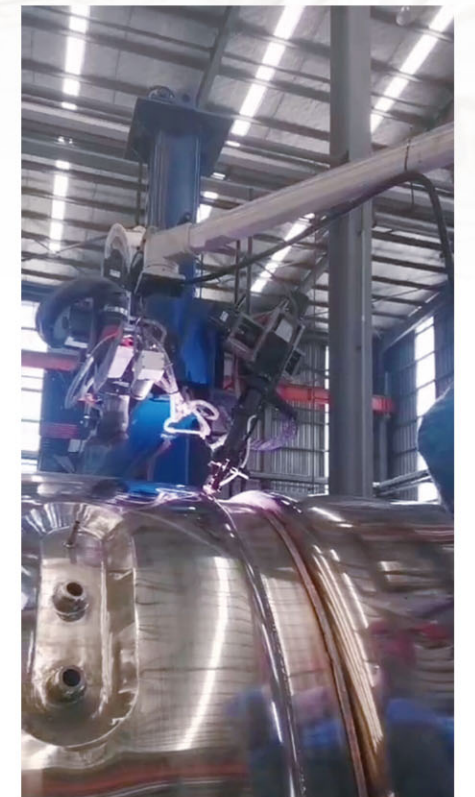Field manufacturing: fermenters with diameter over  $\Phi$ 4,200mm and maximum volume of 1,200m<sup>3</sup>; storage tanks with maximum volume of 10,000m<sup>3</sup>.

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
In biopharmaceutical industry, requirement on fermentation system is more stringent :

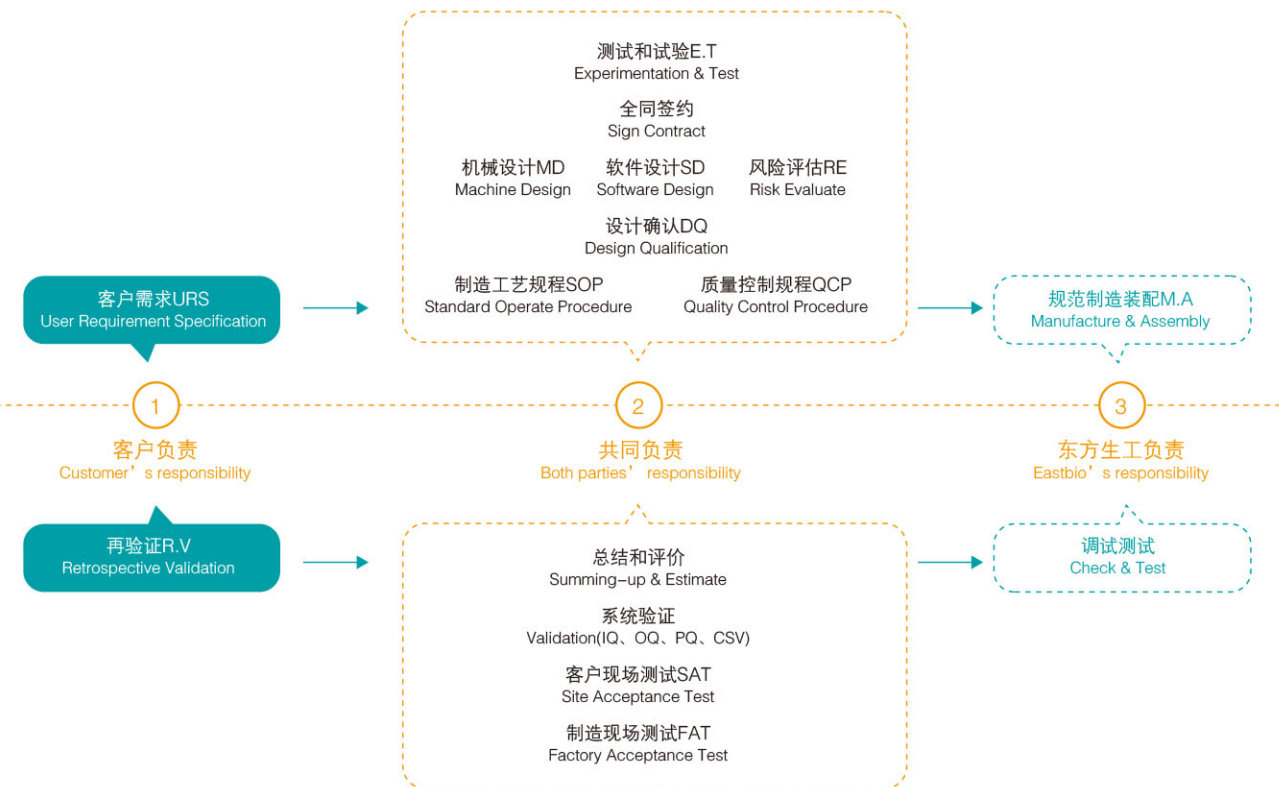All basic features of the equipment must be consistent with GMP and FDA requirements. In terms of equipment for high-end application, Eastbio adopts the international common process in the whole production process to make it to meet the GMP requirements of biomedicine.

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Our professional and efficient after-sales service team provides customers with fault consultation, remote diagnosis and on-site solutions, makes irregular phone calls to understand the operation of equipment, prompt service response while users need it, and always show the service concept of Efficient and Honest.

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EastBio established Jiangsu Multifunctional Fermentation Equipment Engineering Research Center – a professional research and service organization, under the supports of Jiangsu Science and Technology Department, the People's Government of Zhenjiang and Zhenjiang Science and Technology Bureau, which covers the entire process of biotechnological industrialization including strain optimization, fermentation, extraction, filling and packing, etc.

Jiangsu Multifunctional Fermentation Equipment Engineering Research Center has invested RMB 60 million Yuan in building a pilot base, which has a 7000 square meters' construction area and comprises of following facilities: Strain Cultivation Area: designed according to the U.S. FDA standard, there are 4 sterile rooms, 3 culturing rooms, 1 inspection room, and 3 strain preservation rooms;

QC Area: designed according to the U.S. FDA standard, there are 3 sterile laboratories (checking total bacteria counts, disease bacteria, fungi and yeast) and 1 endotoxin room – equipped with imported advanced physical and chemical analysis equipment such as ultra violet analyzer, high-performance liquid chromatography, gas chromatographic analyzer, infrared spectrometer, atomic absorption spectrometer and so on.

Complete Fermentation Laboratory: with 4 sets of 5L automatic fermenters, 3 sets of 20L-30L two levels automatic fermenters, 1 set of 50L-500L fully automatic pilot fermenters.

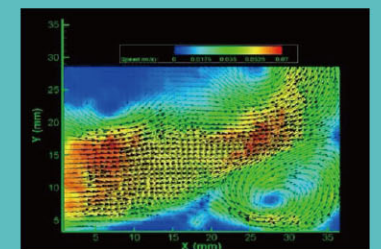
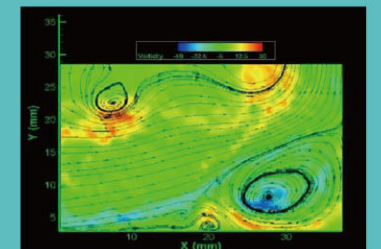
Automatic Fermentation Pilot Plant: equipped with 2 sets of 500L seed fermenter, 2 sets of 3000L fermenters, 2 sets of 5000L fermenters, 2 sets of 20KL fermenters and supporting medium preparation and feeding tanks.

Downstream Extraction Equipment: complete sets of equipment for high-speed centrifugal separation, filter press, solvents, hydrolysis, precipitation and drying, etc.

Ancillary Utility System: solvent recovery, RO water, wastewater treatment, nitrogen extraction, compressed air and steam etc.

Process Automation Laboratory: R&D of bioprocess automation, which involves 13 process parameters, including CER, OUR and so on.

Fermentation Agitation System Research Laboratory: including design of agitation system, simulation of fluid CFD, 3-D printing of agitation system, checking and test of agitation system, and balance test of agitation system.

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
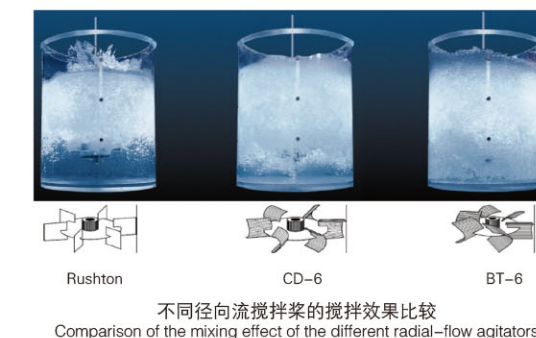
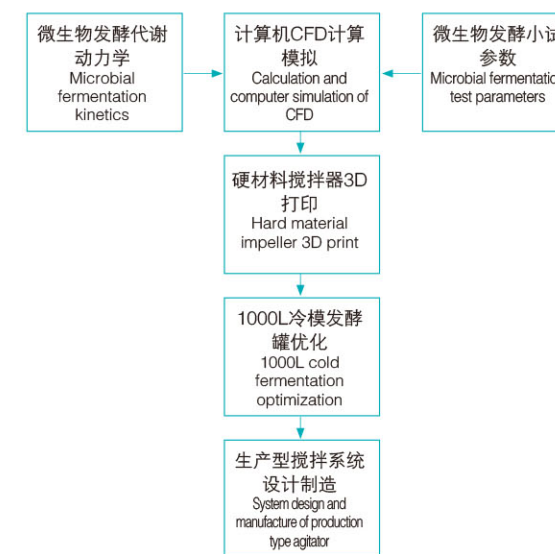
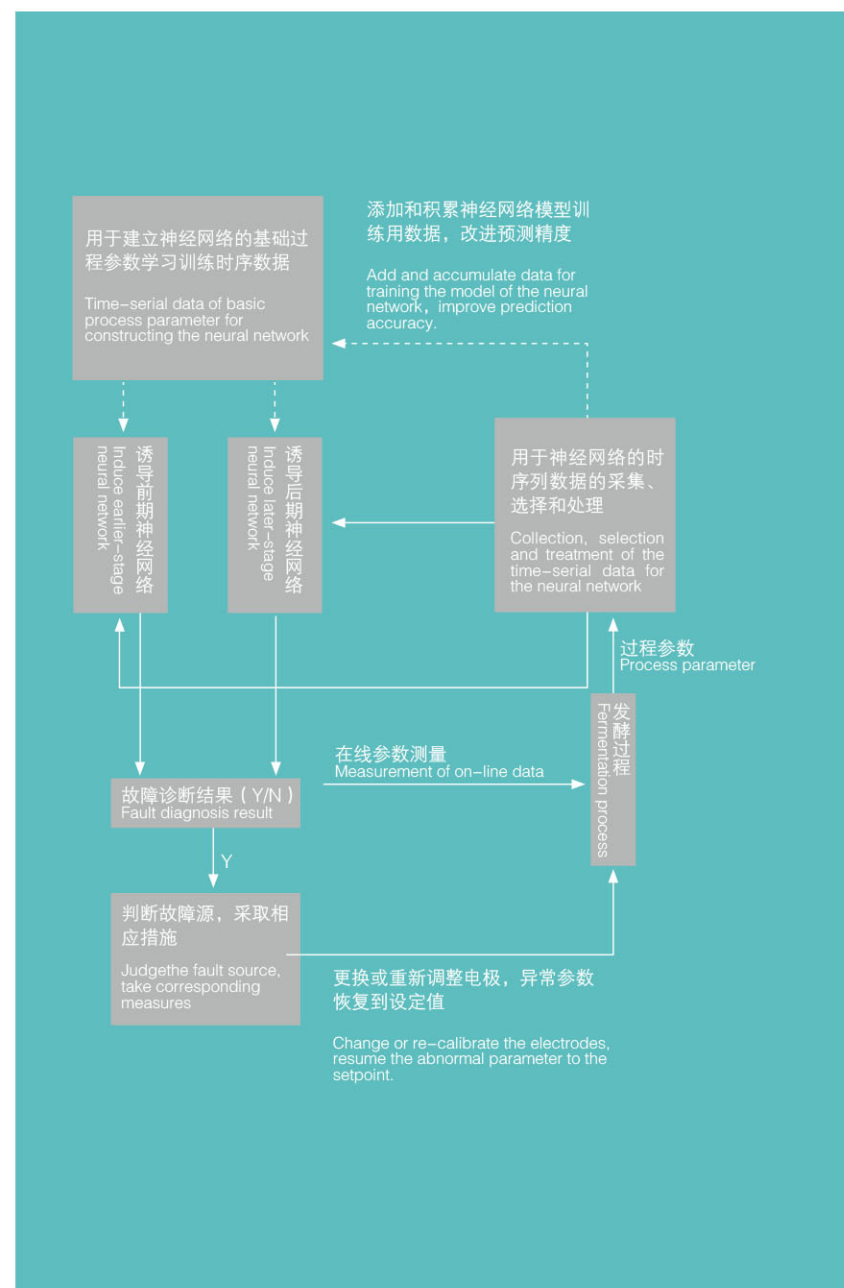
### Intelligent Automation Technology in Bioprocess

As a life body, microorganism shows, in the process of growth and metabolism, dynamic and complicated characteristics, i.e. polymorphic phase, multi-scale, spatial and temporal transient. These characteristics make it very difficult for offline or simple PID control system to adapt to the real-time changes of physiological state of the microorganisms in fermentation process. Artificial intelligence-based automatic control system is a fuzzy-logic control mode. It has played an increasingly important role in the theory and practice of modern fermentation. Eastbio constructed adaptive-control modules and a fault-diagnosis expert system to develop a fully automatic control system and software kits, by adopting method of online detection of process parameters, analysis technology of secondary parameters and artificial intelligent control theory, and integrating process monitoring of physiological status and online-analysis technology of metabolic flux. Now, the control system of Eastbio is able to detect and control more than 13 parameters.

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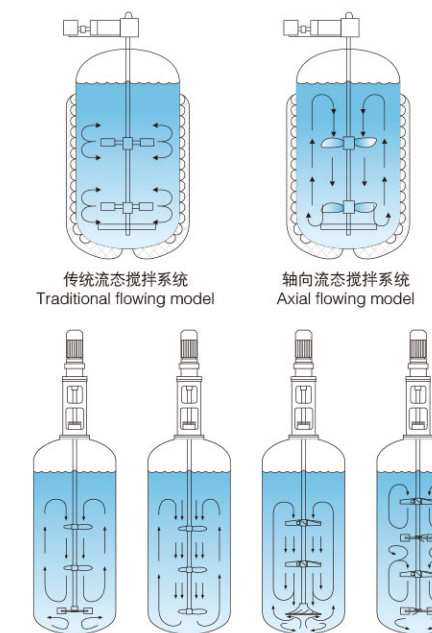
- JE: 轴向流节能搅拌系统  
Energy-saving axial flow agitation system
- JA: 高密度培养大补料工艺搅拌系统  
Mixing system for mass feeding process of high-density culture
- JS: 适合丝状微生物, 粗醪发酵  
Agitation system for filamentous microorganism culture and fermentation with solid material
- JT: 高溶氧搅拌系统  
Agitation system with high dissolved oxygen

### Mass Transfer and Mixing Technology in Bioprocess

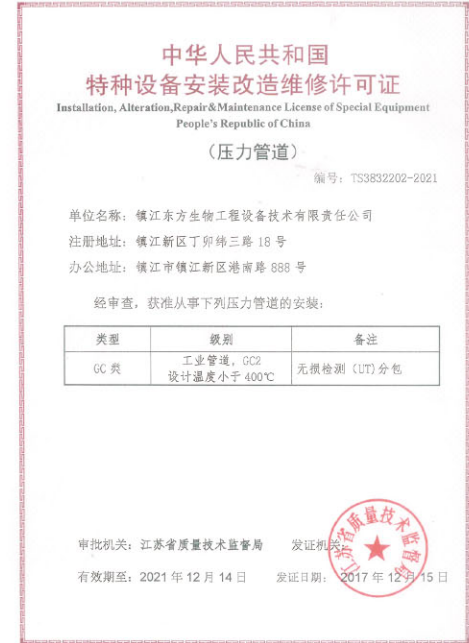Effective mixing and mass transfer technology is crucial for large-scale aerobic fermentation. It is an important factor that directly affects respiratory metabolism of microorganisms and energy-saving benefit of fermentation process. On the basis of traditional scaling-up experience, Eastbio explores options to apply computational technology of fluid dynamics in the design of large-scale aerobic fermentation process. The agitation system and the structure of fermenter can be optimized by mimicking with computer the fluid-flow field, shear-stress field, gas-holdup distribution, energy dissipation, distribution of volumetric mass transfer coefficient and biological reaction changing process in the fermenter. Then the principle and method for rational scaling-up in fermentation process could be established

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METTLER TOLEDO Company Authorization  
(In China, Eastbio uses various sensors and instruments from METTLER TOLEDO, most renowned sensor making company in the world)

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