Operation Manual


Component wR

Large Capacity Centrifuge

Product Model Name:	
---------------------	--

Date of Purchase:



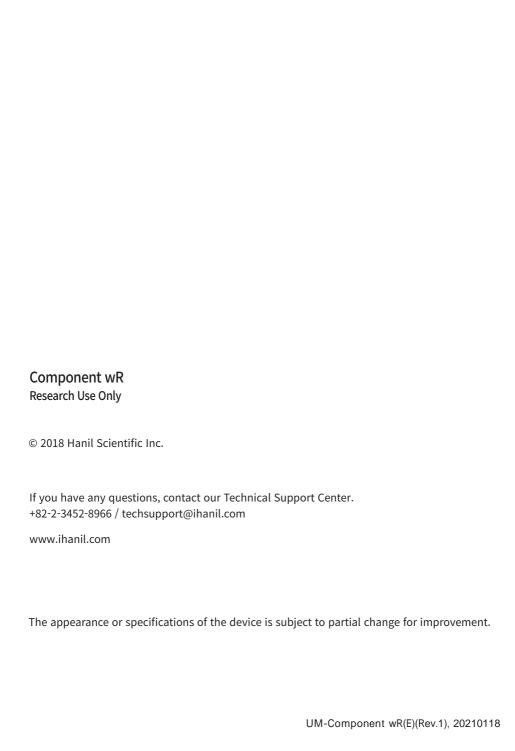

INDEX

1. General Considerations	4
1.2 Transport & Storage - 5	
1.3 Safety label on the Device - 5	
1.4 Electric safety - 5	
2. Product Description	6
2.1 Structure - 6	
2.2 Delivery package - 6	
2.3 Technical Specifications - 7	
3. Installation	8
3.1 Packing Inspection- 8	
3.2 Installation - 8	
3.3 Power Connection - 10	
3.4 Opening/Closing the Lid - 10	
3.5 Using the Rotor - 11	
3.6 Loading Tubes - 12	
3.7 Recommended Tubes - 13	
4. Operation	14
4.1 Control Panel - 14	17
4.2 Setting Speed - 15	
4.3 Setting Run Time - 16	
4.4 Setting Temperature - 16	
4.5 Setting Temperature Limit - 16	
4.6 Setting Acceleration/Deceleration - 17	
4.7 Saving/Calling up a Program - 17	
4.8 Start/Stop/Pulse Run - 18	
4.0 Start/Stop/Fulse Rull 10	
5. Maintenance	19
5.1 Care Instructions - 19	
5.2 Cleaning - 19	
5.3 Care Instructions for Refrigerated device - 20	
6. Troubleshooting	21
8	21
6.1 Possible Problems - 21	
6.2 Error Codes - 22	
6.3 Emergency Lid Open - 23	
7. Rotor	24

1. General Considerations

1.1 Safety

Follow precautions and all the safety requirements described on this user manual to prevent any damage and failure of device and loss of lives.

- 1. The device should be installed on flat surface to maintain level.
- 2. Check the voltage to be used, before connecting the device to the power source.
- Only use rotors, parts, and accessories provided by Hanil Scientific Inc.
 Hanil Scientific Inc. is not responsible for damages of the device and accidents caused by using parts and accessories not recommended.
- 4. Do not exceed the maximum rated speed of the rotor or buckets in use.
- 5. Make sure to prepare necessary safety measures before using samples that are toxic or radioactive samples or pathogenic or samples or infectious blood.
- 6. Substances that may generate volatile or explosive vapor can not be deviced.
- 7. The balancing work of samples should done in advance before operation.
- 8. To ensure safe use of the device, do not expose the device to strong acids, strong bases, cesium, salt, or alkaline detergents.
- 8. If the device is contaminated by toxic or radioactive samples or infectious blood samples, remove contaminants completely and take needful actions such as ventilation or isolation of device.
- 9. Before operation, rotor and chamber should be dry.
- 10. Do not attempt to slow or stop the spinning rotor by hand.
- 11. Only device with rotor and rotor door firmly tightened.
- 12. Do not block vents.
- 13. When serving the device, be sure to remove contaminants in advance.
- 14. Please contact the place of purchase or Hanil Scientific Inc. for product repairs.
- 15. According to IEC61010-2-020 maintain a 30cm "clearance envelope" around the device while the rotor is spinning.
- 16. Turn the power switch off after using the device.
- 17. Unplug the power plug before cleaning or left unused for a long period of time.
- 18. Don't lean against the device

1. General Considerations

1.2 Transport & Storage



- The device and the accessories must only be stored in dry rooms.
- Only lift and transport the device with sufficient number of helpers.

-Storage-

Ambient temperature 5°C~40°C Maximum relative humidity 30%~85% Air pressure 500~1060hpa

-Transport-

Ambient temperature -10°C~40°C Maximum relative humidity 10%~90% Air pressure 500~1060hpa

1.3 Safety label on the Device



Insert the rotor and tubes symmetrically.

Firmly tighten the rotor door.

Watch your fingers when close the lid.



Mark indicating danger and warning.



Mark indicating a place in danger of electric shock.




Mark indicating earth grounding

1.4 Electric safety

- 1. It is recommended that switchgear or circuit breaker and overcurrent protection devices should be installed near the device.
- 2. Use a power cord only provided with the device.
- 3. Do not use an extension cord.
- 4. Do not place anything on the powercable.
- 5. If you have the following emergencies, shut off the power supply and unplug the powercord from outlet and contact your place of purchase or Hanil Scientific Inc.
- Unusual noises or smell from the equipment.
- Damage or wear of a power cord.
- Breakdown of circuit breaker, fuse or safety device.
- If you spill liquid on the equipment.
- If the equipment has been damaged.

2. Product Description

2.1 Structure

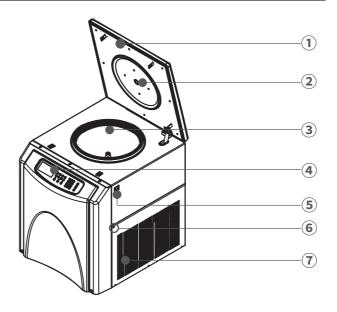

- 1) Lid
- ② RPM measuring window
- ③ Chamber
- 4 Control Panel
- **5** Power Switch
- **6** Emergency Release Hole
- 7 Ventilation Slots

2.2 Delivery package

- Main Device
- Operating manual
- T-wrench
- Rotors & Accessories on request

2. Product Description

2.3 Technical Specifications

Max. RPM (Swing-out Rotor)	4,500 rpm
Max. RCF (Swing-out Rotor)	5,096 xg
Time	< 100 hr, continuous, pulse
Max. Capacity (Swing-out Rotor)	6 x 750 mL
Temperature range	-20°C to 40°C
ACC/DEC steps	10/10 steps
Program memory	100
Rotor Identification	Automatic
Imbalance cutoff / tracking	Yes
Noise level	< 65 dB
Dimension (W x D x H, mm)	668.7 x 800.6 x 891.3
Weight without rotor	230 kg
Power requirement	3.0 kW
Power input (V, Hz)	210~240V 60 Hz (110V, 50 Hz optional)
Cat. No.	CP-wR

3.1 Packing Inspection



- · Check packing conditions carefully, before unpacking.
- Contact Hanil Scientific Inc. immediately if damages found.
- Check the delivery for completeness.

▶ You can get contact details on the packing box or manual.

3.2 Installation

3.2.1 Selecting the Location

Installation on hard and flat ground.

- Centrifuge should be installed on hard and flat place.
- If the centrifuge is installed in an inclined place, the shaft may be bent due to the weight of the rotor.

Good ventailation.

• For air circulation and safety, maintain a 30cm "clearance envelope" around the centrifuge while the rotor is spinning.



Constant temperature/humidity

- Centrifuge equipped with the sensitive electronic software which is fragile with humidity and temperature.
- Must avoid direct ray or heater and be put in the ambience of controlled temp.and air.

Avoid the corrosive gas

- Install the centrifuge in a place where corrosive gas is not generated.
- Sulfur dioxide gas and chlorine gas may cause corrosion.

Leveling

• The device should be operated on a flat surface. Use a leveler tool to check the device is level.

3.2 Installation


3.2.1 Leveling the Device



- · If the device is not installed horizontally, it may cause vibration, noise or malfunction.
- 1. Place the device on a hard and flate place.
- 2. Put the leveler tool on the device.



- 2. Turn the leveling foot on the botton of the device to level the device.
- ▶ Adjust the leveling foot so that the bubble in the leveler tool is at the center of the line.

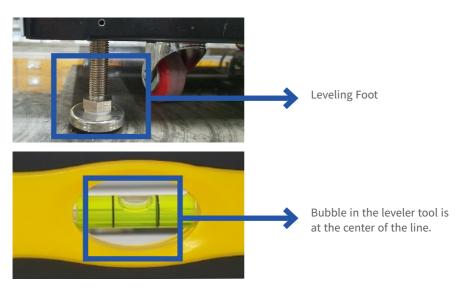

Technical Support: +82-2-3452-8966 / techsupport@ihanil.com | 9/24

3.3 Power Connection



- Connect the device to voltage sources which correspond to the electrical requirements on the label attached to the device.
- Only use a socket proviced Hanil Scientific Inc.
- 1. You can choose a embedded socket or a movable socket
- 2. Switch the centrifuge on using the power switch key on the right side of the device.

3.4 Opening/Closing the Lid



• If the lid is not closed, the device does not operate.

Opening the lid

1. Press the Lid button

Closing the lid

1. Put both hands on the lid and close it with light force.

3.5 Using the Rotor



• Do not disassemble the rotor by user.

This device dose not necessary to remove the rotor after use. Do not disassemble the rotor by the user.

3.6 Loading Tubes



· Use recommended tubes by Hanil Scientific Inc.

- · Always use the same type of tube.
- · Weight, Density and volume of opposing tubes must be same.
- · Check symmetric loading by balancing the tubes with scales.







3.7 Recommended Tubes

Swing-out Rotor: WS750-6


Adaptor	Bore (Ø x L, mm)	Tube	Tube Capacity	Tube Material	Max. g-force(xg)	Tube dimension O.D. x h, mm
-	99 x 103 (B750)	Thermo 75003699	750 mL	Polypropylene	-	98 x 133
TR500(750)	75.5 x 98.7	Hanil HI_TB-w500	500 mL	PP		73.7 x 143
TR500c(750)	99 x 58	Corning 431123	500 mL	PP	6,000 xg	95.5 x 148.6
TR250(750)	62.3 x 87	Nalgene 3120-0250 Nalgene 3122-0250 Autofil 1801-RLS Autofil1803-RLS Autofil 1802-RLS Autofil 1804-RLS	250 mL conical	PPCO PC PP PP PC PC	13,200 xg 27,500xg 8,250xg 27,500xg 8,250xg 31,700xg	61.8 x 125.2 61.8 x 125.2 61 x 127 61 x 127 61 x 127 61 x 127
TR50-7	29.2 x 97	Nalgene 3118-0050 Nalgene 3119-0050	50 mL	PC PPCO	50,000xg 100,605xg	28.8 x 106.4 28.8 x 105.9
TR50c-5	29 x 91	FALCON 352070	50 mL conical	PP	16,000xg	30 x 115
		BD vacutainer 7.5mL~10mL Φ16mm x h 75~100mm	7.5 ~10 mL	-	-	16 x 75~100
TR15-19	17.2 x 87	vacuette 7~9 mL tubes Φ 16mm x h 100mm	7~ 9 mL	-	-	16 x 100
		Nalgene 3110-0150 (cap available separately DS3111-0016)	15 mL	PPCO	50,000 xg	15.8 x 112.8
T15c-14	17 x 89	SPL 50015 Eppendorf 0030122151	15 mL conical	PP PP	9,000 xg 19,500 xg	17 x 120 17 x 121
T10 01	16 07	BD vacutainer 7.5~10 mL Φ16mm x h 75~100mm	7.5 ~10 mL	-	-	16 x 75~100
T10-21 16 x 87	vacuette 7~9 mL tubes Φ 16mm x h 100mm	7~ 9 mL	-	-	16 x 100	
TE 24	13.2 x 60 (3 mL)	BD 1.8ml~7 mL vacutainer Φ13mm x h 75~100mm	1.8~7 mL	-	-	13 x 75~100
	13.2 x 87 (5 mL)	vacuette 1~ 6 mL tubes Φ13mm x h 75~100mm	1~ 6 mL	-	-	13 x 75~100

Tube Material

PPCO: Polypropylene copolymer

PP: Polypropylene PC: Polycarbonate

4.1 Control Panel

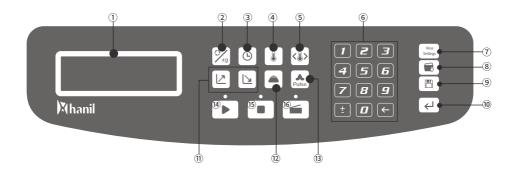

	Button	Description
1)	Display	- Indicate Speed/Time/Temperature/Acc.Dec.
2	Speed	- Press once: Setting the RPM / Press twice: Setting the RCF
3	Time	- Setting run time
4	Temp	- Setting temperature
(5)	Temp Limit	- Setting temperature limit range
6	Numeric buttons	- Enter a value using the numeric buttons
7	View Setting	- Verifying the setting parameters during centrifugation
8	Program call	- Saving a program
9	Program save	- Calling up a program
10	Enter	- Press the enter after setting the parameters
11)	Acceleration (↗) Deceleration (↘)	- Setting acceleration step - Setting deceleration step
(12)	Rotor ID.	- When rotor id is not identified automatically, enter the rotor ID. manually
13)	Pulse	- Short spin
14)	Start	- Starting centrifugation
(15)	Stop	- Stopping centrifugation
16)	Lid	- Openg the lid

4.2 Setting Speed

[Setting RPM]

- 1. Press the Speed button once.
 - ▶ RPM setting mode: RPM(rpm) display flashes.
- 2. Press the numeric buttons to set the desired RPM value.
 - ▶ Up to 4,500rpm
- 3. Press the Enter button.

[Setting RCF]

- 1. Press the Speed button twice.
 - ► RCF setting mode: RCF(xg) display flashes.
- 2. Press the numeric buttons to set the desired RCF value.
 - ▶ Up to 5,096xg
- 3. Press the Enter button.

4.3 Setting Run Time

- 1. Press the Time button.
 - ▶ Time setting mode: Time 00:00:00 display flashes.
- 2. Press the numeric buttons to set the desired run time.
- ▶ Up to <100hr
- 3. Press the Setup/Enter button.

4.4 Setting Temperature

- 1. Press the Temp button.
- 2. Press the numeric buttons to set the desired temperature.
- ► -20°C ~ 40 °C
- 3. Press the Setup/Enter button.

4.5 Setting Temperature Limit

Temperature range that stops operation when the temperature is lower or higher than the set temperature.

- 1. Press the Temp limit button.
- 2. Press the numeric buttons to set the desired temperature limit range.
 - ▶ Default value: ±30°C
- 3. Press the Setup/Enter button.

4.6 Setting Acceleration/Deceleration

- 1. Press the ACC or DEC button.
 - ► Acceleration setting mode: ACC (↗) display flashes.
 - ▶ Deceleration setting mode: DEC (\(\sigma\)) display flashes.
- 2. Press the numeric buttons to set the desired ACC or DEC rate.
- ► ACC rate: 0~9 / DEC rate: 0~9
- 3. Press the Enter button.

4.7 Saving/Calling up a Program

[Saving]

- 1. Set the RPM(or RCF) / Run time / Accel, Decel rate / Temperature
- 2. Press the Program Save button.
- 3. Press the numeric buttons to set the desired program number.
 - ▶ Up to 0~99:100
- 4. Press the Enter button.

[Calling up]

- 1. Press the Program Call button.
- 2. Press the numeric buttons to set the desired program number.
- 3. Press the Enter button.

4.8 Start/Stop/Pulse Run

[Start a run]

- 1. Check all parameters are correct.
- 2. Press the Start button.

[Stop a run]

1. Press the Stop button.

[Pulse]

It is for quick and short spin down.

After reaching the set speed, the centrifugation is decelated and stopped.

- 1. Set the desired speed.
 - ► Refer to [4.2 Setting Speed]
- 2. Press the Pulse button.

5. Maintenance

5.1 Care Instructions

- 1. Regularly inspect the rotor chamber for check the motor shaft is normal.
- 2. Regularly check the rotor and accessories to be sure there is no damage.
- 3. Rotate the motor shaft with your hand to make sure it turns smoothly.
- 4. To prevent the rotor from sticking, lubicate the rotor drive hole which contacts motor shaft regularly.

5.2 Cleaning



· Before cleaning the centrifuge, be sure to switch off the device and disconnect the power cord.

[Oustside of the device]

- 1 Clean the outside of deivice with a soft and dry cloth.
- ▶ If the device is contaminated, use a mild cleaning fluid to clean.
- 2. Do not use aggressive chemicals on the device such as alchol, benzene, acetone or phenol.
- 3. Make sure do not scratch the surface of deivice when cleaning it.
 - ▶ Do not use a metal sponge.
- ▶ If the device is rusted, remove it with a mild detergent and wipe it with a dry cloth.

[Chamber]

- 1. If the rotor chamber is not dry, wipe moisture from the chamber with a dry cloth.
- 2. If the rotor chamber is dirty, clean it with a mild fluid to clean.

[Rotor]

- 1. After centrifugation, remove the rotor from the chamber and keep it upside down.
- 2. If any sample is spilt inside the rotor, clean it and dry well.

[Disposal]

1. In case of product is to be disposed of, the local wastes laws and regulations are to be observed.

5. Maintenance

5.3 Care Instructions for Refrigerated device

[Open the lid after centrifugation]


- 1. Wipe up the moisture in the chamber with a soft cloth.
- 2. Leave the centrifuge lid open after centrifugation.

[Remove the condensation water]

1. Remove the drain cap and empty the condensation water regularly.

[Remove dust]

1. Remove any dust from the ventilation slots regularly.

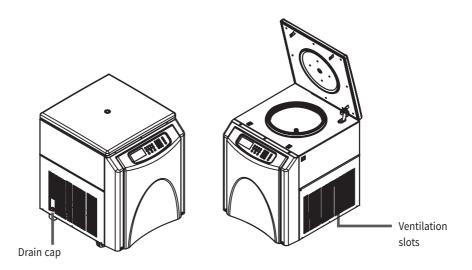

6. Troubleshooting

6.1 Possible Problems

Problems	Recommended Action		
Power failure	Check the power cord connection.Check the power fuse of the device.		
Centrifugation cannot be started	- Check the lid is closed completely.		
Lid cannot be opened	- Press the 'Lid open button'. - Check the power connection.		
Lid cannot be closed	- Remove the dirt at the door latch and close the lid Check the lid latch is not damaged.		
	- Check the device is installed on the hard and flat place		
Unusual noise and vibration	- Reinstall the rotor symmetrically Reinstall the tubes symmetrically.		
Imbalance error message is displayed	7000rpm 0:30:0 28°1 Imbalance Error! - If there is a problem with the imbalance sensor, an imbalance error message is displayed In this case, pressing the start button does not work Contact Hanil techsupport.		

6. Troubleshooting

6.2 Error Codes


Error	Message	Cause	Recommended Action
E1	Imbalance error	•Imbalance is detected.	 Check the tightness of rotor loading. Check the proper loading of tubes. Check the stable ground or worktable.
E2	Over speed error	10% over the set RPM.Incorrect tuning of motor and controller.	Check controller and motor. Restart the device
E3	Motor start error	• Temp. rises inside of the motor.	Restart the device.
E4	Door Open! Fast stop!	- When the lid is opened during operation.	Restart the device
E5	Low Speed error	• Can not reach 100rpm within 2sec. after start the run.	Restart the device

6. Troubleshooting

6.3 Emergency Lid Open

If the device lid cannot be opened, you can activte the emergency open manually.

- 1. Wait for rotor to stop before activating the emergency open.
- 2. Insert the T-wrench into the emergency open hole and press it until the lid is opened.

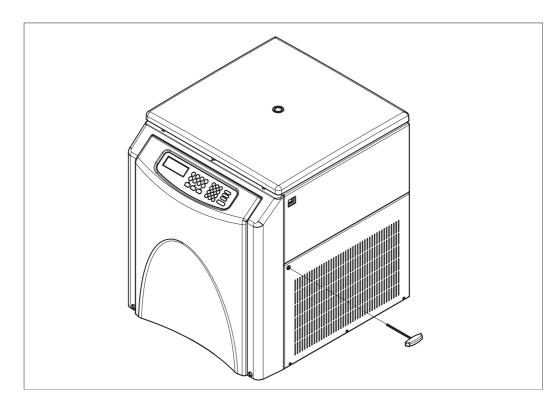

7. Rotors and Accessories

WS750-6





Round Bucket B750 B750S (Cap available)











∠90° Max. RPM: 4,500 Size (W x D x H)

: Ø262 x 262 mm (w/o windshield)









Adaptor TR5-24

Bucket	Required Adaptor	Tube Capacity Bottom Type	Tube per Adaptor / Rotor	Bore (Ø x L, mm)	Max. height for tube fit (mm) (w/ cap)	Max. height for tube fit (mm) (w/o cap)	Max. RPM (rpm) Max. RCF (xg)
	-	750 mL Flat	1/4	99 x 103	129	161.5	4,500 5,096
	TR500(750)	500 mL Flat	1/4	75.5 x 98.7	125.2	129	4,500 5,042
	TR500c(750)	500mL conical Conical	1/4	99 x 58	126.7	154.3	4,500 5,096
	TR250(750)	250 mL Flat	1/4	62.3 x 87	124.5	149	4,500 5,003
B750S /	TR50-7	50 mL Round	7 / 28	29.2 x 97	124	131.8	4,500 5,003
B750	TR50c-5	50 mL conical Conical	5 / 20	29 x 91	116.2	125.5	4,500 5,096
	TR15-19	15 mL Round	19 / 76	17.2 x 87	114.6	122.4	4,500 5,003
	TR15c-14	15 mL conical Conical	14 / 56	17 x 89	117	124.8	4,000 3,479
	TR10-21	10 mL Round	21 / 84	21 / 84	113.9	121.7	4,500 5,003
	TR5-24	3 mL / 5 mL Round	24 / 96	13.2 x 60(3 mL) 13.2 x 87(5 mL)	113.7	121.5	4,500 5,003
TM96(750-4))	MTP	1/4	88 x 128.5 x 42			4,500 4,345



hanil

16, Arayuk-ro, Gimpo, REP of Korea +82-2-3452-8966 technicalsupport@ihanil.com www.ihanil.com